



# Quick Guide Omegon® Microstar Camera Installation

## 1. Installing the Omegon® Microstar Camera Driver


Click on ->Omegon Microstar Camera Setup.exe file from the CD-ROM file

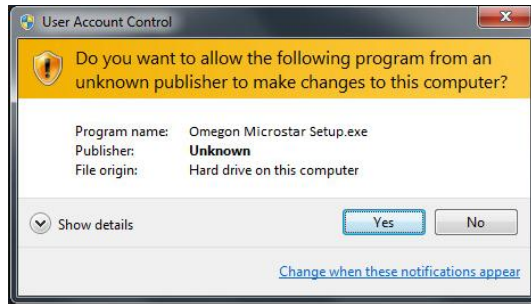


The User Account Control message will show up. Click 'Yes' to continue.

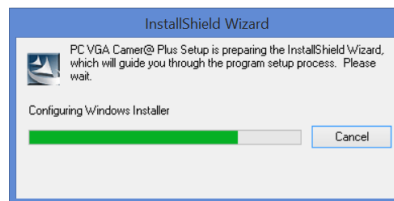

The PC VGA Camer@ plus Setup Driver will start to install

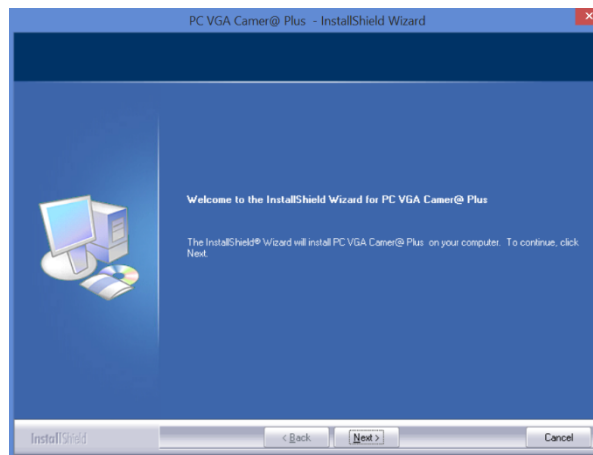

Click ->Next to start

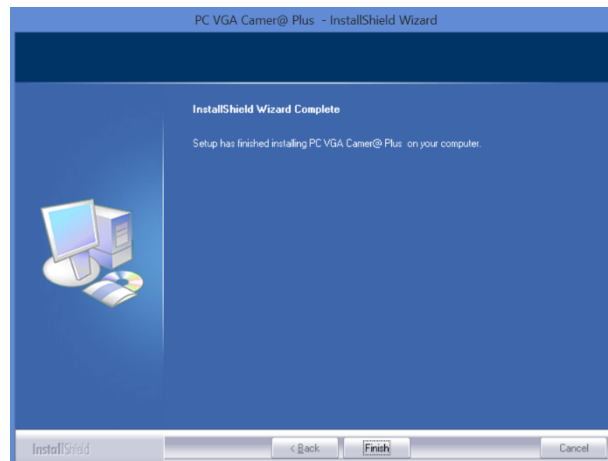


Click -> Finish when install is complete


## CONNECT THE CAMERA USB CABLE TO THE COMPUTER

### 2. Using the supplied capture video Omegon® AMcap.exe


Click on ->Omegon® AMcap.exe file from the CD-ROM file

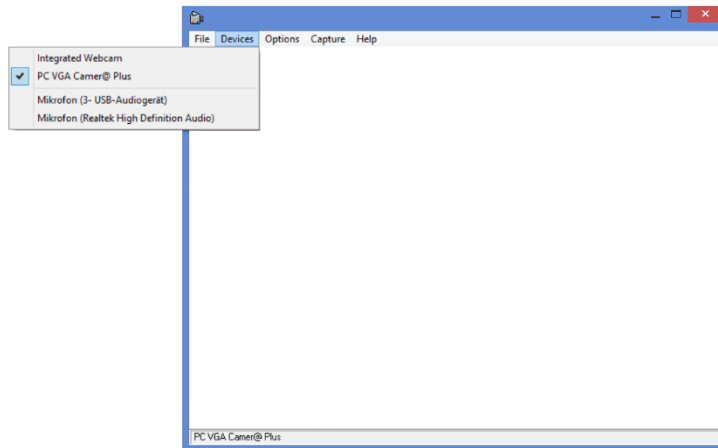


Select the Microstar camera from the available camera list. **Devices->Options>PC VGA Camer@ Plus**

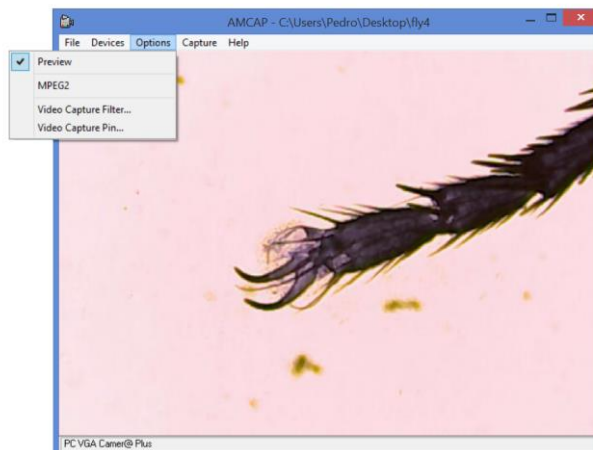


Select **Options->Preview**.

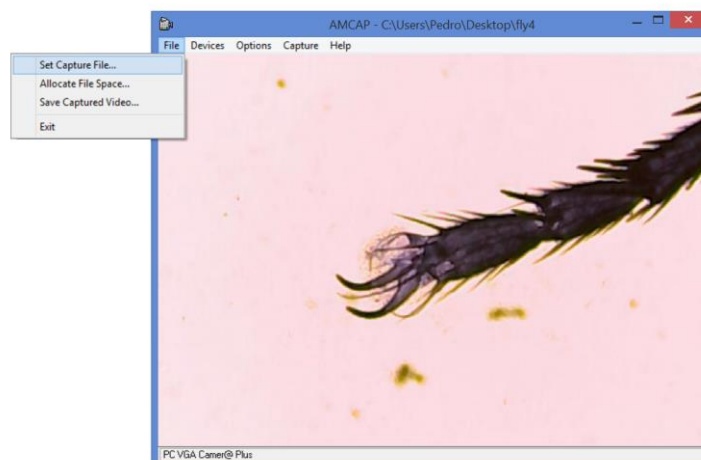


To save a capture select the File name. Click **File-> Set Capture File**  
Choose a name for the file (notice: the file is an .AVI file if the extension is not assigned to the file please do it manually. The file show be of the type File.AVI)

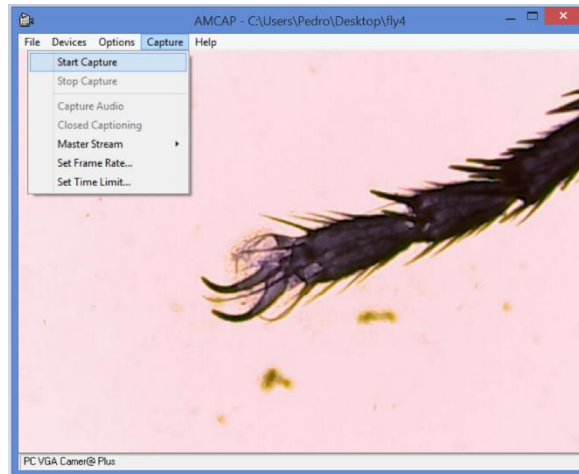

Click **Capture->Start Capture**

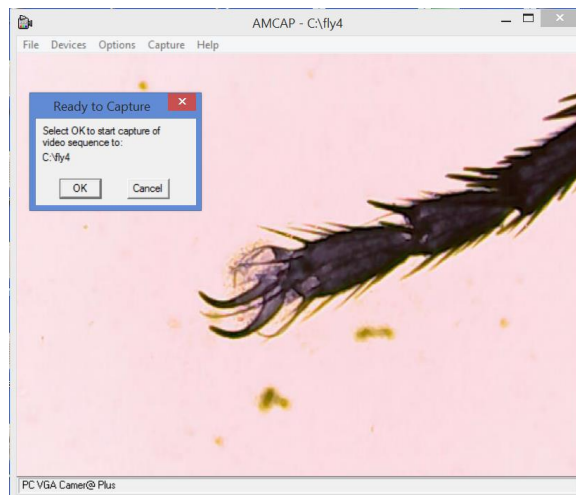

Check if file path is correct and press '**OK**' to start capture.

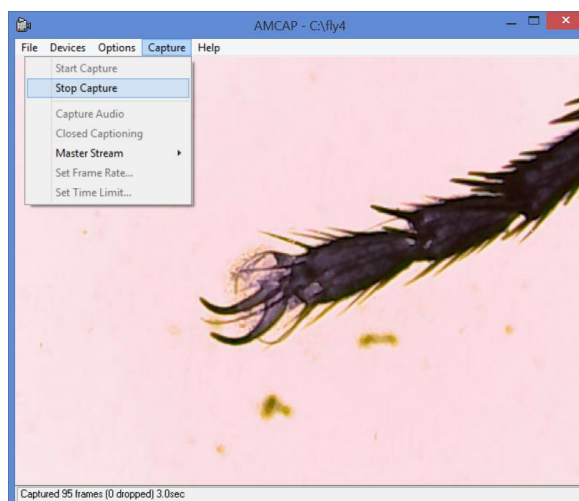

Click **Capture->Stop Capture** to end capture.